

## Swivel-Mounted Multi-Function-Head - ONE-TWIST


This multi-function-head will be found only with our automatic and semi-automatic flash-point testers.

Which are:

**ABA 4** (acc. Abel Method)

**TAG 4** (acc. TAG Method)

**PMA 4, PMP 4 and PM 4** (acc. Pensky-Martens Method).

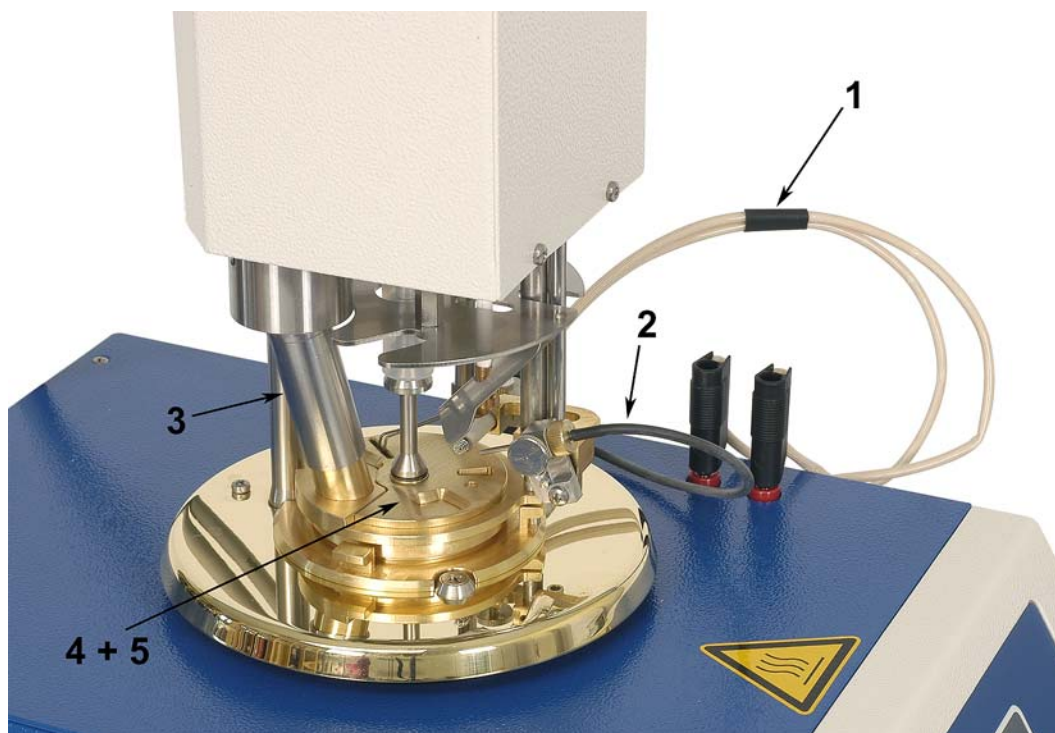

### Description of the Multi-Function-Head

Figure with gas-igniter.

The gas-ignition will be lit up by the electric-igniter automatically.

1. Electric-Igniter
2. Gas-Igniter
3. Multi-Detector
4. Cup
5. Cover

**Unbelievable fast swivel-around process within 2 steps only:**



**Step 1:**

Simply insert the complete test insert (cup with cover and multi-detector).

Please use the tongs (supplied with the instrument) or the optional handle which will be attached to the cup permanently.


**Step 2:**

**Snap-in the head - Ready!:**

Just lift up the head a little bit and turn it clockwise.

All connections with stirrer, cup opening, igniter position, temperature- & flash-point indicator are made simultaneously after twisting the head into position.

It's so simple - only with us!

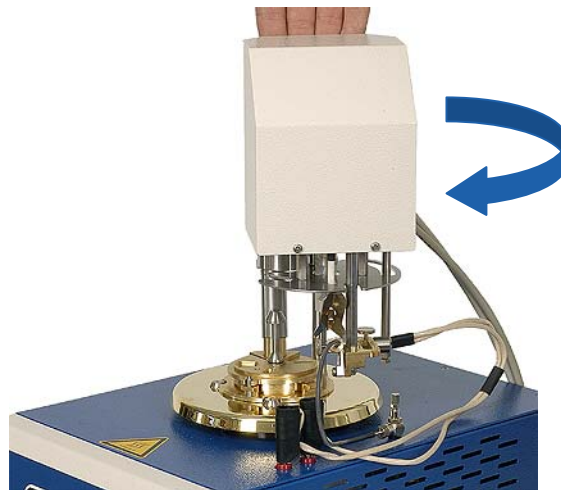

The next Flash-Point Test can be started.

Figure with electric-ignition.  
The procedure is the same with gas-ignition.

**Compare this with competing instruments**